



CERTIFICATION by CUSTODIAN of RECORDS

Tucson Police Department Crime Laboratory
1306 W. Miracle Mile, Tucson, AZ 85705

I hereby certify that these documents are a true and correct copy of the periodic maintenance* for the Intoxilyzer 8000 Serial Number 80-000323 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

TSy
11-3-2015

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 10/15/2015 Cal 1 0.101 AC Cal 2 0.101 AC

Between the dates of 9/21/2015 and 10/15/2015 the instrument was operating properly and accurately.

The instrument was taken out of service for repair on
The instrument was returned to service on

William L Loomis

William L. Loomis


10/22/2015

ADPS QAS Permit # 33720

SUBSCRIBED AND SWORN to before me this 22nd, day of October, 2015.

Elma Escarcega

Notary Public

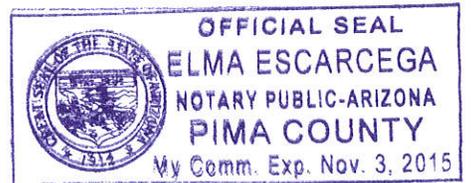

*Includes documents specified in A.A.C. R13-10-104.A.5

EXHIBIT G-4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

STANDARD QUALITY ASSURANCE PROCEDURE

1 of 9
WLL
TSh
11-3-2015

QA SPECIALIST W. LOOMIS 53545 AGENCY Tucson Police Department

DATE 10/15/2015 TIME 16:17

INTOXILYZER SERIAL # 80-000323 LOCATION CRIME LAB

(✓) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- (✓) 1. Clock time check.
- (✓) 2. Date check.

OPERATIONAL TESTS

- (✓) 1. Alcohol-free subject test result 0. 000 AC.
- (✓) 2. Error recognition logic system functioning.
Not a Successfully Completed Test Sequence printed
- (✓) 3. Proper sample recognition system.
Not a Successfully Completed Test Sequence printed
Deficient sample printed.
- (✓) 4. Standard Calibration Check standard 0. 100 AC
Result: 0. 101 AC

Instrument is operating properly and accurately. YES NO

COMMENTS

SIGNATURE Will Loomis

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-000323
Core Version: 8105.48
10/15/2015

16:20:46

Standard Lot#: 12415100A4
Last Changed By: W. LOOMIS #53545

QAS: W. LOOMIS #53545
TUCSON PD

Operator: W. LOOMIS #53545
TUCSON PD

Subject: SQAP, TEST
DOB: 11/11/1987
Sex: M Weight: 222

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	16:22:03
Diagnostic Test	Pass	16:22:37
Air Blank	0.000	16:23:04
0.100 Cal Check	0.101	16:23:25
Air Blank	0.000	16:23:54
Subject Test	0.000	16:24:12
Air Blank	0.000	16:24:39
Wait	ABT*	
Air Blank	0.000	16:25:17

*Sequence Aborted

2 of 9
WLL
Bh

Not a Successfully Completed Test Sequence

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-000323
Core Version: 8105.48
10/15/2015

16:25:52

3 of 9
WLL
TBY

Standard Lot#: 12415100A4
Last Changed By: W. LOOMIS #53545

QAS: W. LOOMIS #53545
TUCSON PD

Operator: W. LOOMIS #53545
TUCSON PD

Subject: SQAP, TEST
DOB: 11/11/1987
Sex: M Weight: 222

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	16:28:18
Diagnostic Test	Pass	16:28:52
Air Blank	0.000	16:29:19
0.100 Cal Check	RFI*	16:29:33
Air Blank	IPS**	16:29:41

*RFI Detect
**Improper Sample

Not a Successfully Completed Test Sequence

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-000323
Core Version: 8105.48
10/15/2015

16:30:51

Standard Lot#: 12415100A4
Last Changed By: W. LOOMIS #53545

QAS: W. LOOMIS #53545
TUCSON PD

Operator: W. LOOMIS #53545
TUCSON PD

Subject: SQAP, TEST
DOB: 11/11/1987
Sex: M Weight: 222

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	16:32:23
Diagnostic Test	Pass	16:32:57
Air Blank	0.000	16:33:24
0.100 Cal Check	0.102	16:33:45
Air Blank	0.000	16:34:15
Subject Test	0.255*	16:37:22
Air Blank	0.000	16:37:55
Air Blank	0.000	16:38:23
Subject Test	XXX**	16:38:44
Air Blank	0.000	16:39:18

*Deficient Sample

**Invalid Sample

Begin new deprivation period

4 of 9
WLL
BY

Not a Successfully Completed Test Sequence

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-000323
Core Version: 8105.48
10/15/2015 16:40:11

Standard Lot#: 12415100A4
Last Changed By: W. LOOMIS #53545

QAS: W. LOOMIS #53545
TUCSON PD

Operator: W. LOOMIS #53545
TUCSON PD

Test	g/210L	Time
Air Blank	0.000	16:40:41
0.100 Cal Check	0.101	16:41:02
Air Blank	0.000	16:41:31

Preventative Maintenance
Performed -- Timer Reset

5 of 9
WLL
Bh

EXHIBIT G-2

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

6 of 9
WLL
By
11-3-2015

QA Specialist W. LOOMIS 53545 AGENCY Tucson Police Department
DATE 10/15/2015 TIME 16:17
Intoxilyzer Serial # 80-000323 Location CRIME LAB

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.
OR
Pour a standard alcohol concentration solution _____ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C.
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 101 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Will Loomis

DPS Form Exh G-2 (Rev. 05-01)

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-000323
Core Version: 8105.48
10/15/2015 16:17:56

7 of 9
WLL
BH

Standard Lot#: 12415100A4
Last Changed By: W. LOOMIS #53545

QAS: W. LOOMIS #53545
TUCSON PD

Operator: W. LOOMIS #53545
TUCSON PD

Test	g/210L	Time
Air Blank	0.000	16:18:26
0.100 Cal Check	0.101	16:18:46
Air Blank	0.000	16:19:16

Preventative Maintenance
Performed -- Timer Reset



TUCSON POLICE DEPARTMENT
CRIME LABORATORY

INTOXILYZER SERVICE RECORD
FOR
CALIBRATION GAS TANK CHANGE

8 of 9
WLL
By
11-3-2015

Date: 10/1/2015
 QAS: W. LOOMIS 53545
 Intoxilyzer Serial Number: 80-000323
 New Lot Number: 12415100 A4
 New Expiration Date: 6/5/2017

- (✓) 1. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 2. Remove the old gas tank and install the new one.
- (✓) 3. Hit the "ESC" key twice and swipe the QAS card.
- (✓) 4. Go to the "M-Maintenance" menu and select "S-Standard Change." Select "D" for dry gas and enter the following information:
 Enter Target Value: 0.100
 Enter Control Lot #: From the tank label
 Enter Exp Date: From the tank label (mm/dd/yyyy)
 Edit info Y/N? Y or N
 Update Std Info? Y ENTER
 "Please wait saving..."
- (✓) 5. Hit the "ESC" key twice.
- (✓) 6. Go to the "M-Maintenance" menu and select "T-Control Testing." Select "D" for dry control test. After selection is made press ENTER.
- (✓) 7. Air blank completed.
- (✓) 8. Calibration check completed. Test results 0. 103 AC.
- (✓) 9. Air blank completed.
- (✓) 10. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Will Loomis

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-000323
Core Version: 8105.48
10/01/2015 15:53:23

Standard Lot#: 12415100A4
Last Changed By: W. LOOMIS #53545

QAS: W. LOOMIS #53545
TUCSON PD

Operator: W. LOOMIS #53545
TUCSON PD

Test	g/210L	Time
Air Blank	0.000	15:53:52
0.100 Cal Check	0.103	15:54:13
Air Blank	0.000	15:54:42

Preventative Maintenance
Performed -- Timer Reset

9 of 9
WLL
TS4