



CERTIFICATION by CUSTODIAN of RECORDS  
Tucson Police Department Crime Laboratory

*Handwritten initials*

I hereby certify that these documents are a true and correct copy of the periodic maintenance\* for the Intoxilyzer 8000 Serial Number 80-001165 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 12.1.2006 Cal 1 0.097 AC Cal 2 0.098 AC

The instrument was placed into service on 12.1.2006.


*Terry S. Gallegos*

Terry S. Gallegos

1.4.2007

ADPS QAS Permit # 33169

SUBSCRIBED AND SWORN to before me this 4, day of January, 2007.

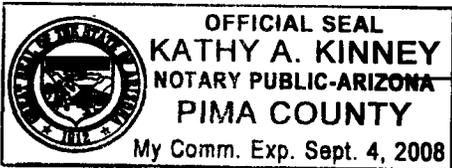

*Kathy A. Kinney*

Notary Public

\*Includes documents specified in A.A.C. R13-10-104.A.5

1058  
D4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD QUALITY ASSURANCE PROCEDURE

QA SPECIALIST Terry Gallegos 28047 AGENCY Tucson Police Department

DATE 12-1-2006 TIME 152

INTOXILYZER SERIAL # 80-001165 LOCATION Lab

( ✓ ) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- ( ✓ ) 1. Clock time check.
- ( ✓ ) 2. Date check.

OPERATIONAL TESTS

- ( ✓ ) 1. Alcohol-free subject test result 0. 000 AC.
- ( ✓ ) 2. Error recognition logic system functioning.  
Not a Successfully Completed Test Sequence printed ✓✓
- ( ✓ ) 3. Proper sample recognition system.  
Not a Successfully Completed Test Sequence printed ✓  
Deficient sample printed. ✓
- ( ✓ ) 4. Standard Calibration Check standard 0. 100 AC  
Result: 0. 098 AC

Instrument is operating properly and accurately. YES  NO

COMMENTS

---



---



---



---

SIGNATURE 

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

2 of 8  
T84

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-001165  
12/01/2006 11:55:44

Standard Lot#: 502602I  
Last Changed By: T. GALLEGOS #28847

QAS: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

Operator: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

Subject: SQAP, TEST  
DOB: 11/11/1111  
Sex: M Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	11:57:41
Diagnostic Test	Pass	11:58:09
Air Blank	0.000	11:58:36
0.100 Cal Check	0.097	11:58:55
Air Blank	0.000	11:59:24
Subject Test	0.000	11:59:51
Air Blank	0.000	12:00:18
Five Minute Wait	ABT*	
Air Blank	0.000	12:00:53

\*Sequence Aborted

**Not a Successfully Completed Test Sequence**

3 of 8  
TG

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT *je*  
Serial Number : 80-001165  
12/01/2006 12:01:06

---

Standard Lot#: 502602I  
Last Changed By: T. GALLEGOS #28847

---

QAS: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

---

Operator: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

---

Subject: SQAP, TEST  
DOB: 11/11/1111  
Sex: M Weight: 111

---

15 Minute Deprivation Period? Yes

---

Test	g/210L	Time
Air Blank	IPS*	12:01:47
Air Blank	0.000	12:02:15

\*Improper Sample

**Not a Successfully Completed Test Sequence**

4 of 8  
TS6

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT *js*  
Serial Number : 80-001165  
12/01/2006 12:02:34

---

Standard Lot#: 502602I  
Last Changed By: T. GALLEGOS #28847

---

QAS: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

---

Operator: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

---

Subject: SQAP,TEST  
DOB: 11/11/1111  
Sex: M Weight: 111

---

15 Minute Deprivation Period? Yes

---

Test	g/210L	Time
Air Blank	0.000	12:03:27
Diagnostic Test	Pass	12:03:55
Air Blank	0.000	12:04:22
0.100 Cal Check	0.099	12:04:41
Air Blank	0.000	12:05:10
Subject Test	RFI*	12:06:51
Air Blank	0.000	12:07:19

\*RFI Detect

**Not a Successfully Completed Test Sequence**

5 of 8  
T4

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-001165  
12/01/2006 12:08:57

Standard Lot#: 502602I  
Last Changed By: T. GALLEGOS #28847

QAS: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

Operator: T. GALLEGOS #28847  
TUCSON PD CRIME LAB

Subject: SQAP,TEST  
DOB: 11/11/1111  
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	12:09:58
Diagnostic Test	Pass	12:10:25
Air Blank	0.000	12:10:53
0.100 Cal Check	0.099	12:11:11
Air Blank	0.000	12:11:41
Subject Test	0.021*	12:14:50
Air Blank	0.000	12:15:20
Air Blank	0.000	12:15:47
Subject Test	XXX**	12:16:34
Air Blank	0.000	12:17:08

\*Deficient Sample  
\*\*Invalid Sample  
Begin new deprivation period

**Not a Successfully Completed Test Sequence**

*Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.*

INTOXILYZER 8000

Location: TUCSON POLICE DEPT

Serial Number: 80-001165

12/01/2006

12:17:40

6078  
TS

-----  
Standard Lot#: 502602I

Last Changed By: T. GALLEGOS #28847

-----  
QAS: T. GALLEGOS #28847

TUCSON PD CRIME LAB  
-----

Test	g/210L	Time
Air Blank	0.000	12:18:07
Cal Check	0.098	12:18:26
Air Blank	0.000	12:18:55

  
-----  
Operator's Signature

EXHIBIT G-2

7 of 8  
TJG

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

jk

STANDARD CALIBRATION CHECK PROCEDURE

QA Specialist Terry Gallegos<sup>28947</sup> AGENCY Tucson Police Department  
DATE 12-1-2006 TIME 1152  
Intoxilyzer Serial # 80-001165 Location Lab

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0.100 AC.  
OR  
Pour a standard alcohol concentration solution \_\_\_\_\_ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0.097 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Terry Gallegos

DPS Form Exh G-2 (Rev. 05-01)

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.


8 of 8  
D4

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number: 80-001165  
12/01/2006 11:52:11

-----  
Standard Lot#: 502602I  
Last Changed By: T. GALLEGOS #28847  
-----

QAS: T. GALLEGOS #28847  
TUCSON PD CRIME LAB  
-----

Test	g/210L	Time
Air Blank	0.000	11:52:38
Cal Check	0.097	11:52:57
Air Blank	0.000	11:53:27

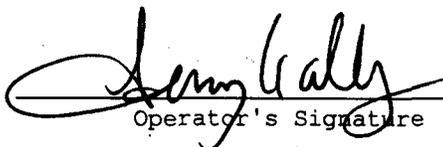  
Operator's Signature



TUCSON POLICE DEPARTMENT  
CRIME LABORATORY

*je*

INTOXILYZER SERVICE RECORD

*1af1  
Bh*

Intoxilyzer Serial Number: 80-001165

Initial Service Date: 12-1-2006

CMI, Inc. Certificate of Calibration Date: 2.4.2005