



CERTIFICATION by CUSTODIAN of RECORDS
Tucson Police Department Crime Laboratory

TSH
12-13-07

I hereby certify that these documents are a true and correct copy of the periodic maintenance* for the Intoxilyzer 8000 Serial Number 80-002264 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 11-19-07 Cal 1 0.101 AC Cal 2 0.099 AC

Between the dates of 10/19/07 and 11/19/07 the instrument was operating properly and accurately.

The instrument was taken out of service for repair on 11/9/07.
The instrument was returned to service on 11/14/07.


Clare Conley

Clare L. Conley

12-7-07

ADPS QAS Permit # 33314

SUBSCRIBED AND SWORN to before me this 7, day of December, 2007.

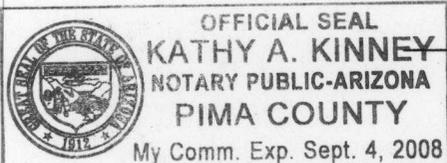

Kathy A Kinney
Notary Public

*Includes documents specified in A.A.C. R13-10-104.A.5

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

EXHIBIT G-4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

1 of 7 cu

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

T84
12.13.07

STANDARD QUALITY ASSURANCE PROCEDURE

QA SPECIALIST Clare Conley AGENCY Tucson Police Department
DATE 11-19-07 TIME 9:36am
INTOXILYZER SERIAL # 80-002264 LOCATION DUI Headquarters

(✓) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- (✓) 1. Clock time check.
- (✓) 2. Date check.

OPERATIONAL TESTS

- (✓) 1. Alcohol-free subject test result 0. 000 AC.
- (✓) 2. Error recognition logic system functioning.
Not a Successfully Completed Test Sequence printed
- (✓) 3. Proper sample recognition system.
Not a Successfully Completed Test Sequence printed
Deficient sample printed.
- (✓) 4. Standard Calibration Check standard 0. 100 AC
Result: 0. 099 AC

Instrument is operating properly and accurately. YES NO

COMMENTS None.

SIGNATURE Clare Conley

INTOXILYZER 8000

Location: TUCSON POLICE DEPT

Serial Number : 80-002264

11/19/2007

09:47:49

2 of 7 cu
134

Standard Lot#: 719801I

Last Changed By: S. SULLIVAN

#34971

QAS: C. CONLEY

#34517

TUCSON PD

Operator: C. CONLEY

#34517

TUCSON PD

Subject: TEST, SQAP

DOB: 11/11/1111

Sex: M

Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	09:48:48
Diagnostic Test	Pass	09:49:15
Air Blank	0.000	09:49:42
0.100 Cal Check	0.101	09:50:01
Air Blank	0.000	09:50:30
Subject Test	0.000	09:50:47
Air Blank	ABT*	09:50:54
Air Blank	0.000	09:51:22

*Sequence Aborted

Not a Successfully Completed Test Sequence

Pursuant to section 17-322 Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number : 80-002264
11/19/2007

09:51:27

387 cu
754

Standard Lot#: 719801I
Last Changed By: S. SULLIVAN #34971

QAS: C. CONLEY #34517
TUCSON PD

Operator: C. CONLEY #34517
TUCSON PD

Subject: TEST, SQAP
DOB: 11/11/1111
Sex: M Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	RFI*	09:52:05
Air Blank	IPS**	09:52:24

*RFI Detect

**Improper Sample

Not a Successfully Completed Test Sequence

Pursuant to section 28-3204 Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number : 80-002264
11/19/2007 09:53:08

457cm
TS4

Standard Lot#: 719801I
Last Changed By: S. SULLIVAN #34971

QAS: C. CONLEY #34517
TUCSON PD

Operator: C. CONLEY #34517
TUCSON PD

Subject: TEST, SQAP
DOB: 11/11/1111
Sex: M Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	09:53:59
Diagnostic Test	Pass	09:54:27
Air Blank	0.000	09:54:54
0.100 Cal Check	0.101	09:55:13
Air Blank	0.000	09:55:42
Subject Test	0.297*	09:58:51
Air Blank	0.000	09:59:23
Air Blank	0.000	09:59:51
Subject Test	XXX**	10:00:06
Air Blank	0.000	10:00:40

*Deficient Sample

**Invalid Sample

Begin new deprivation period

Not a Successfully Completed Test Sequence

Pursuant to section 322A of the Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000
Location: TUCSON POLICE DEPT
Serial Number: 80-002264
11/19/2007 10:01:07

5067 cu
TS4

Standard Lot#: 719801I
Last Changed By: S. SULLIVAN #34971

QAS: C. CONLEY #34517
TUCSON PD

Test	g/210L	Time
Air Blank	0.000	10:01:35
Cal Check	0.099	10:01:53
Air Blank	0.000	10:02:23

Clare Conley

Operator's Signature

EXHIBIT G-2

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY
STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000
STANDARD CALIBRATION CHECK PROCEDURE

687 cu
By
12.13.07

QA Specialist Clare Conley AGENCY Tucson Police Department
DATE 11-19-07 TIME 9:36 am
Intoxilyzer Serial # 80-002264 Location DUI Headquarters

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.
OR
Pour a standard alcohol concentration solution _____ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 101 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Clare Conley

DPS Form Exh G-2 (Rev. 05-01)

INTOXILYZER 8000

Location: TUCSON POLICE DEPT

Serial Number: 80-002264

11/19/2007

09:36:47

7 of 7 cu

Standard Lot#: 719801I

Last Changed By: S. SULLIVAN #34971

134

QAS: C. CONLEY #34517

TUCSON PD

Test g/210L Time

Air Blank 0.000 09:37:14
Cal Check 0.101 09:37:33
Air Blank 0.000 09:38:02

Clare Conley

Operator's Signature



TUCSON POLICE DEPARTMENT
CRIME LABORATORY

INTOXILYZER SERVICE RECORD

Dy Bg
12.13.07 10f2

Date: 11/14/2007

QAS: Terry Gallegos 28847

Intoxilyzer Serial Number: 80-002264

Reason For Service: Breath hose not heating. Officer Hawkins indicated the breath hose problem on 11-9-07.

Date Out of Service: N/A

Calibration Check completed prior to removal from service

Date: N/A

SQAP completed prior to removal from service

Date: N/A

Corrective Action: Sent to CMI, Inc.

Replaced breath hose and returned to service on 11.14.2007.

INTOXILYZER 8000

Location: TUCSON POLICE DEPT

Serial Number: 80-002264

11/14/2007

13:48:12

2 of 2
B4 B4

Standard Lot#: 719801I

Last Changed By: S. SULLIVAN #34971

QAS: S. CIONE #50616

TUCSON PD

Test	g/210L	Time
Air Blank	0.000	13:48:39
Cal Check	0.100	13:48:58
Air Blank	0.000	13:49:27

Sarah Cione

Operator's Signature