



CERTIFICATION by CUSTODIAN of RECORDS

Tucson Police Department Crime Laboratory  
1306 W. Miracle Mile, Tucson, AZ 85705

I hereby certify that these documents are a true and correct copy of the periodic maintenance\* for the Intoxilyzer 8000 Serial Number 80-002759 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

sc  
11/19/13

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 10.30.2013 Cal 1 0.097 AC Cal 2 0.097 AC

Between the dates of — and — the instrument was operating properly and accurately.

The instrument was taken out of service for repair on —.  
The instrument was returned to service on 10.30.2013.

Terry Gallegos

Terry S. Gallegos


11.7.2013

ADPS QAS Permit # 33169

SUBSCRIBED AND SWORN to before me this 7, day of November, 2013.

A. M. Hernandez

Notary Public

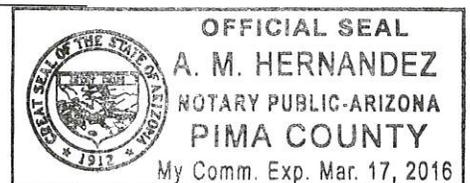

\*Includes documents specified in A.A.C. R13-10-104.A.5



TUCSON POLICE DEPARTMENT  
CRIME LABORATORY

INTOXILYZER SERVICE RECORD

SC  
11/19/13

QAS: Terry Gallegos 28847

Intoxilyzer Serial Number: 80-002759

Date Returned to Service: 10-30-2013

EXHIBIT G-4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD QUALITY ASSURANCE PROCEDURE

10 of 9  
By SC  
11/19/13

QA SPECIALIST Terry Gallegos 28847 AGENCY Tucson Police Department

DATE 10.30.2013 TIME 1738

INTOXILYZER SERIAL # 80-002759 LOCATION CrimeLab

( ✓ ) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- ( ✓ ) 1. Clock time check.
- ( ✓ ) 2. Date check.

OPERATIONAL TESTS

- ( ✓ ) 1. Alcohol-free subject test result 0. 000 AC.
- ( ✓ ) 2. Error recognition logic system functioning.  
Not a Successfully Completed Test Sequence printed
- ( ✓ ) 3. Proper sample recognition system.  
Not a Successfully Completed Test Sequence printed  
Deficient sample printed.
- ( ✓ ) 4. Standard Calibration Check standard 0. 100 AC  
Result: 0. 097 AC

Instrument is operating properly and accurately. YES  NO

COMMENTS

---



---



---



---

SIGNATURE 

INTOXILYZER 8000  
Location: TUCSON PD  
Serial Number: 80-002759  
Core Version: 8105.48  
10/30/2013

17:41:52

2059  
TS4 SC

---

Standard Lot#: 30412100A1  
Last Changed By: T. GALLEGOS #28847

---

QAS: T. GALLEGOS #28847  
TUCSON PD

---

Operator: T. GALLEGOS #28847  
TUCSON PD

---

Subject: TEST,SQAP  
DOB: 11/11/1987  
Sex: F Weight: 111

---

15 Minute Deprivation Period? Yes

---

Test	g/210L	Time
Air Blank	0.000	17:42:49
Diagnostic Test	Pass	17:43:24
Air Blank	0.000	17:43:52
0.100 Cal Check	0.098	17:44:13
Air Blank	0.000	17:44:42
Subject Test	0.000	17:45:02
Air Blank	0.000	17:45:29
Wait	ABT*	
Air Blank	0.000	17:46:03

\*Sequence Aborted

**Not a Successfully Completed Test Sequence**

INTOXILYZER 8000  
Location: TUCSON PD  
Serial Number: 80-002759  
Core Version: 8105.48  
10/30/2013

17:46:10

30f 9  
TS4  
5c

---

Standard Lot#: 30412100A1  
Last Changed By: T. GALLEGOS #28847

---

QAS: T. GALLEGOS #28847  
TUCSON PD

---

Operator: T. GALLEGOS #28847  
TUCSON PD

---

Subject: TEST,SQAP  
DOB: 11/11/1987  
Sex: F Weight: 111

---

15 Minute Deprivation Period? Yes

---

Test	g/210L	Time
Air Blank	0.000	17:47:00
Diagnostic Test	Pass	17:47:35
Air Blank	0.000	17:48:02
0.100 Cal Check	RFI*	17:48:16
Air Blank	IPS**	17:48:29

\*RFI Detect

\*\*Improper Sample

**Not a Successfully Completed Test Sequence**

INTOXILYZER 8000  
Location: TUCSON PD  
Serial Number: 80-002759  
Core Version: 8105.48  
10/30/2013

17:48:37

4 of 9  
B4  
5c

---

Standard Lot#: 30412100A1  
Last Changed By: T. GALLEGOS #28847

---

QAS: T. GALLEGOS #28847  
TUCSON PD

---

Operator: T. GALLEGOS #28847  
TUCSON PD

---

Subject: TEST, SQAP  
DOB: 11/11/1987  
Sex: F Weight: 111

---

15 Minute Deprivation Period? Yes

---

Test	g/210L	Time
Air Blank	0.000	17:49:32
Diagnostic Test	Pass	17:50:07
Air Blank	0.000	17:50:35
0.100 Cal Check	0.098	17:50:56
Air Blank	0.000	17:51:25
Subject Test	0.028*	17:54:33
Air Blank	0.000	17:55:05
Air Blank	0.000	17:55:32
Subject Test	XXX**	17:55:58
Air Blank	0.000	17:56:32

\*Deficient Sample

\*\*Invalid Sample

Begin new deprivation period

**Not a Successfully Completed Test Sequence**

INTOXILYZER 8000  
Location: TUCSON PD  
Serial Number: 80-002759  
Core Version: 8105.48  
10/30/2013 17:57:07

5089  
TS4  
5C

-----  
Standard Lot#: 30412100A1  
Last Changed By: T. GALLEGOS #28847  
-----

QAS: T. GALLEGOS #28847  
TUCSON PD  
-----

Operator: T. GALLEGOS #28847  
TUCSON PD  
-----

Test	g/210L	Time
Air Blank	0.000	17:57:37
0.100 Cal Check	0.097	17:57:58
Air Blank	0.000	17:58:27

Preventative Maintenance  
Performed -- Timer Reset

EXHIBIT G-2

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

6 of 9  
TS4  
SC  
11/14/13

QA Specialist Terry Gallegos 28847 AGENCY Tucson Police Department  
DATE 10-30-2013 TIME 1738  
Intoxilyzer Serial # 80-002759 Location Crime Lab

- ( ✓ ) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.  
OR  
Pour a standard alcohol concentration solution \_\_\_\_\_ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- ( ✓ ) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- ( ✓ ) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- ( ✓ ) 4. Air blank completed.
- ( ✓ ) 5. Calibration check completed. Test results 0. 097 AC.
- ( ✓ ) 6. Air blank completed.
- ( ✓ ) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE 

DPS Form Exh G-2 (Rev. 05-01)

INTOXILYZER 8000  
Location: TUCSON PD  
Serial Number: 80-002759  
Core Version: 8105.48  
10/30/2013 17:38:07

7 of 9  
TGA  
5<

-----  
Standard Lot#: 30412100A1  
Last Changed By: T. GALLEGOS #28847  
-----

QAS: T. GALLEGOS #28847  
TUCSON PD  
-----

Operator: T. GALLEGOS #28847  
TUCSON PD  
-----

Test	g/210L	Time
Air Blank	0.000	17:38:45
0.100 Cal Check	0.097	17:39:06
Air Blank	0.000	17:39:35

Preventative Maintenance  
Performed -- Timer Reset



TUCSON POLICE DEPARTMENT  
CRIME LABORATORY

INTOXILYZER SERVICE RECORD  
FOR  
CALIBRATION GAS TANK CHANGE

80F9  
T54  
SC  
11/19/13

Date: 10.30.2013  
QAS: Terry Gallegos 28847  
Intoxilyzer Serial Number: 80-002759  
New Lot Number: 30412100A1  
New Expiration Date: 12.1.2014

- (✓) 1. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 2. Remove the old gas tank and install the new one.
- (✓) 3. Hit the "ESC" key twice and swipe the QAS card.
- (✓) 4. Go to the "M-Maintenance" menu and select "S-Standard Change."  
Select "D" for dry gas and enter the following information:  
  - Enter Target Value: 0.100
  - Enter Control Lot #: From the tank label
  - Enter Exp Date: From the tank label (mm/dd/yyyy)
  - Edit info Y/N? Y or N
  - Update Std Info? Y ENTER
  - "Please wait saving..."
- (✓) 5. Hit the "ESC" key twice.
- (✓) 6. Go to the "M-Maintenance" menu and select "T-Control Testing."  
Select "D" for dry control test. After selection is made press ENTER.
- (✓) 7. Air blank completed.
- (✓) 8. Calibration check completed. Test results 0. 097 AC.
- (✓) 9. Air blank completed.
- (✓) 10. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Terry Gallegos

INTOXILYZER 8000  
Location: TUCSON PD  
Serial Number: 80-002759  
Core Version: 8105.48  
10/30/2013 17:32:50

9 of 9  
TS4  
SC

-----  
Standard Lot#: 30412100A1  
Last Changed By: T. GALLEGOS #28847  
-----

QAS: #

-----  
Operator: T. GALLEGOS #28847  
TUCSON PD CRIME LAB  
-----

Test	g/210L	Time
Air Blank	0.000	17:33:17
0.100 Cal Check	0.097	17:33:38
Air Blank	0.000	17:34:07

# Certificate of Calibration

This is to certify the calibration of **Intoxilyzer**® serial number 80-002759, manufactured by CMI Inc., a subsidiary of MPD, Inc. of Owensboro, Kentucky, was tested and found to conform to the National Highway Traffic Safety Administration (NHTSA) Standard for Devices to Measure Breath Alcohol (Federal Register, Vol.58, No.179, pp 48705-48710, Sept. 17, 1993) for accuracy and precision. Reference materials are traceable through the National Institute of Standards and Technology (NIST) to the International System of Units (SI).

Date July 12, 2013 Signed Dale Rehner  
Technician

**CMI** INC.

316 East 9<sup>th</sup> Street  
Owensboro, KY 42303  
USA

Part No. 650517 Rev.A

Pursuant to Section 28.1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.



# Certificate of Calibration

This is to certify the calibration of **Intoxilyzer**® serial number 80-002759, manufactured by CMI Inc., a subsidiary of MPD, Inc. of Owensboro, Kentucky, was tested and found to conform to the National Highway Traffic Safety Administration (NHTSA) Standard for Devices to Measure Breath Alcohol (Federal Register, Vol.58 No.179, pp 48705-48710, Sept. 17, 1993) for accuracy and precision. Reference materials are traceable through the National Institute of Standards and Technology (NIST) to the International System of Units (SI).

Date Oct. 22, 2013 Signed Del Risher

Technician

**EMI** INC.

316 East 9<sup>th</sup> Street  
Owensboro, KY 42303  
USA

Part No. 650517 Rev.A

Pursuant to Section 28.1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

304  
TS4  
SC

