



CERTIFICATION by CUSTODIAN of RECORDS
 Tucson Police Department Crime Laboratory
 1306 W. Miracle Mile, Tucson, AZ 85705

TBJ
 4.14.2014

I hereby certify that these documents are a true and correct copy of the periodic maintenance* for the Intoxilyzer 8000 Serial Number 80-006053 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 3/28/14 Cal 1 0.098 AC Cal 2 0.098 AC

Between the dates of 2/27/14 and 3/28/14 the instrument was operating properly and accurately.

The instrument was taken out of service for repair on -.
 The instrument was returned to service on -.

Sarah Cione

Sarah L. Cione


4/1/14

ADPS QAS Permit # 33355

SUBSCRIBED AND SWORN to before me this 1st, day of April, 2014.

Elma Escarcega

Notary Public

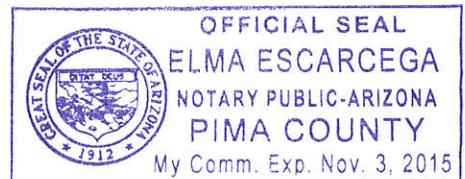

*Includes documents specified in A.A.C. R13-10-104.A.5

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

EXHIBIT G-4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

STANDARD QUALITY ASSURANCE PROCEDURE

1 of 10 sc
TSY
4.18.2014

QA SPECIALIST Sarah Cione AGENCY Tucson Police Department

DATE 3/28/14 TIME 1019

INTOXILYZER SERIAL # 80-006053 LOCATION DUI HQ

(✓) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- (✓) 1. Clock time check.
- (✓) 2. Date check.

OPERATIONAL TESTS

- (✓) 1. Alcohol-free subject test result 0. 000 AC.
- (✓) 2. Error recognition logic system functioning.
Not a Successfully Completed Test Sequence printed
- (✓) 3. Proper sample recognition system.
Not a Successfully Completed Test Sequence printed
Deficient sample printed.
- (✓) 4. Standard Calibration Check standard 0. 100 AC
Result: 0. 098 AC

Instrument is operating properly and accurately. YES NO

COMMENTS SQAP started on 3/27/14 - did cal. check, but then tank pressure not recognized. Performed maintenance, then SQAP performed on 3/28/14.

SIGNATURE Sarah Cione

INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/28/2014

10:21:20

2 of 10_{sc}
B4

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: S. CIONE #50616
TUCSON PD

Operator: S. CIONE #50616
TUCSON PD

Subject: SQAP, SQAP
DOB: 11/11/1980
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 10:22:30 |
| Diagnostic Test | Pass | 10:23:05 |
| Air Blank | 0.000 | 10:23:32 |
| 0.100 Cal Check | 0.099 | 10:23:53 |
| Air Blank | 0.000 | 10:24:23 |
| Subject Test | 0.000 | 10:24:45 |
| Air Blank | 0.000 | 10:25:13 |
| Wait | ABT* | |
| Air Blank | 0.000 | 10:25:52 |

*Sequence Aborted

Not a Successfully Completed Test Sequence

INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/28/2014

10:26:04

3 of 10_{sc}
TSg

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: S. CIONE #50616
TUCSON PD

Operator: S. CIONE #50616
TUCSON PD

Subject: SQAP, SQAP
DOB: 11/11/1980
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 10:27:16 |
| Diagnostic Test | Pass | 10:27:51 |
| Air Blank | 0.000 | 10:28:18 |
| 0.100 Cal Check | 0.099 | 10:28:39 |
| Air Blank | 0.000 | 10:29:09 |
| Subject Test | RFI* | 10:29:39 |
| Air Blank | IPS** | 10:29:48 |

*RFI Detect

**Improper Sample

Not a Successfully Completed Test Sequence

INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/28/2014

10:29:55

4 of 10 SC
TS4

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: S. CIONE #50616
TUCSON PD

Operator: S. CIONE #50616
TUCSON PD

Subject: SQAP, SQAP
DOB: 11/11/1980
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 10:31:08 |
| Diagnostic Test | Pass | 10:31:43 |
| Air Blank | 0.000 | 10:32:10 |
| 0.100 Cal Check | 0.099 | 10:32:31 |
| Air Blank | 0.000 | 10:33:01 |
| Subject Test | 0.310* | 10:36:08 |
| Air Blank | 0.000 | 10:36:41 |
| Air Blank | 0.000 | 10:37:09 |
| Subject Test | XXX** | 10:37:41 |
| Air Blank | 0.000 | 10:38:15 |

*Deficient Sample

**Invalid Sample

Begin new deprivation period

Not a Successfully Completed Test Sequence

INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/28/2014 10:38:30

5 of 10
sc
TS4

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: S. CIONE #50616
TUCSON PD

Operator: S. CIONE #50616
TUCSON PD

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 10:38:58 |
| 0.100 Cal Check | 0.098 | 10:39:19 |
| Air Blank | 0.000 | 10:39:49 |

Preventative Maintenance
Performed -- Timer Reset

EXHIBIT G-2

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

6 of 10
SC
TSY
4.14.2014

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

QA Specialist Sarah Cione AGENCY Tucson Police Department
DATE 3/28/14 TIME 1019
Intoxilyzer Serial # 80-006053 Location DUIHQ

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.
OR
Pour a standard alcohol concentration solution _____ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 098 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Sarah Cione

DPS Form Exh G-2 (Rev. 05-01)

INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/28/2014 10:19:35

7 of 10
SC

TS4

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: S. CIONE #50616
TUCSON PD

Operator: S. CIONE #50616
TUCSON PD

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 10:20:04 |
| 0.100 Cal Check | 0.098 | 10:20:25 |
| Air Blank | 0.000 | 10:20:54 |

Preventative Maintenance
Performed -- Timer Reset

EXHIBIT G-2

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

8 of 10^{sc}
TSN
4.14.2014

QA Specialist Sarah Cione AGENCY Tucson Police Department
DATE 3/27/14 TIME 851
Intoxilyzer Serial # 80-006053 Location DUI HQ

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.
OR
Pour a standard alcohol concentration solution _____ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 099 AC.
- (✓) 6. Air blank completed.
- () 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Sarah Cione

DPS Form Exh G-2 (Rev. 05-01)

INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/27/2014 08:51:01

9 of 10
5
TS4

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: S. CIONE #50616
TUCSON PD

Operator: S. CIONE #50616
TUCSON PD

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 08:51:29 |
| 0.100 Cal Check | 0.099 | 08:51:50 |
| Air Blank | 0.000 | 08:52:20 |

Preventative Maintenance
Performed -- Timer Reset



TUCSON POLICE DEPARTMENT
 CRIME LABORATORY
 INTOXILYZER SERVICE RECORD

10 of 10_{sc}

TS4
 4.14.2014

Date: 3/28/14

QAS: Sarah Cione

Intoxilyzer Serial Number: 80-006053

Reason For Service: Instrument does not recognize any tank pressure, even with full tank. Subject test cannot proceed.

Date Out of Service: N/A

Calibration Check completed prior to removal from service

Date: -

SQAP completed prior to removal from service

Date: -

Corrective Action: Sent to CMI, Inc.

Replaced regulator



TUCSON POLICE DEPARTMENT
CRIME LABORATORY

1 of 2 sc

INTOXILYZER SERVICE RECORD
FOR
CALIBRATION GAS TANK CHANGE

TS4
4.14.2014

Date: MARCH 21st 2014

QAS: A. MARQUEZ #47674

Intoxilyzer Serial Number: 80-006053

New Lot Number: 25913100 A2

New Expiration Date: 10-1-2015

OS
OS
OS

1. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
2. Remove the old gas tank and install the new one.
3. Hit the "ESC" key twice and swipe the QAS card.
4. Go to the "M-Maintenance" menu and select "S-Standard Change." Select "D" for dry gas and enter the following information:

Enter Target Value: 0.100
 Enter Control Lot #: From the tank label
 Enter Exp Date: From the tank label (mm/dd/yyyy)
 Edit info Y/N? Y or N
 Update Std-Info? Y ENTER
 "Please wait saving..."

OS
OS

5. Hit the "ESC" key twice.
6. Go to the "M-Maintenance" menu and select "T-Control Testing." Select "D" for dry control test. After selection is made press ENTER.

OS
OS
OS

7. Air blank completed.
8. Calibration check completed. Test results 0. 099 AC.
9. Air blank completed.

OS

10. Remove and sign printed record. Attach the record to the completed checklist.

SIGNATURE

A. Marquez #47674

2 of 2 sc

TS4



INTOXILYZER 8000
Location: TUCSON PD
Serial Number: 80-006053
Core Version: 8105.48
03/21/2014 17:33:01

Standard Lot#: 25913100A2
Last Changed By: A. MARQUIS #47674

QAS: A. MARQUIS #47674
TUCSON PD

Operator: A. MARQUIS #47674
TUCSON PD

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 17:33:35 |
| 0.100 Cal Check | 0.099 | 17:33:55 |
| Air Blank | 0.000 | 17:34:25 |

MARQUIS 47674

Preventative Maintenance
Performed -- Timer Reset

