



CERTIFICATION by CUSTODIAN of RECORDS

Tucson Police Department Crime Laboratory  
270 S. Stone Avenue Tucson, AZ 85701-1917

TJH  
4.8.2010

I hereby certify that these documents are a true and correct copy of the periodic maintenance\* for the Intoxilyzer 8000 Serial Number 80-000332 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 4-1-10 Cal 1 0.098 AC Cal 2 0.097 AC

Between the dates of 3-4-10 and 4-1-10 the instrument was operating properly and accurately.

The instrument was taken out of service for repair on N/A.  
The instrument was returned to service on N/A.


Clare Conley  
Clare L. Conley

4-1-10

ADPS QAS Permit # 33314

SUBSCRIBED AND SWORN to before me this 1<sup>st</sup> day of April, 2010.

Isabel C. Trujillo  
Notary Public

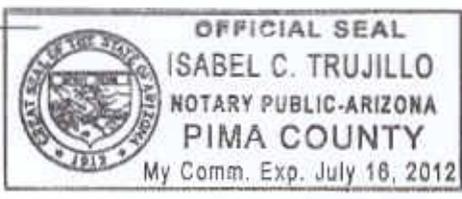

\*Includes documents specified in A.A.C. R13-10-104.A.5

EXHIBIT G-4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD QUALITY ASSURANCE PROCEDURE

1567  
Ch  
TH  
482010

QA SPECIALIST Clare Conley AGENCY Tucson Police Department

DATE 4-1-10 TIME 0837

INTOXILYZER SERIAL # 80-000332 LOCATION ODW DUI HQ

(✓) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- (✓) 1. Clock time check.
- (✓) 2. Date check.

OPERATIONAL TESTS

- (✓) 1. Alcohol-free subject test result 0. 000 AC.
- (✓) 2. Error recognition logic system functioning.  
Not a Successfully Completed Test Sequence printed
- (✓) 3. Proper sample recognition system.  
Not a Successfully Completed Test Sequence printed  
Deficient sample printed.
- (✓) 4. Standard Calibration Check standard 0. 100 AC  
Result: 0. 097 AC

Instrument is operating properly and accurately. YES  NO

COMMENTS None.  
\_\_\_\_\_  
\_\_\_\_\_  
\_\_\_\_\_

SIGNATURE Clare Conley

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-000332  
04/01/2010 08:38:37

2567

OK

BH

Standard Lot#: 642845  
Last Changed By: C. CONLEY #34517

QAS: C. CONLEY #34517  
TUCSON PD

Operator: C. CONLEY #34517  
TUCSON PD

Subject: TEST, SQAP  
DOB: 11/11/1111  
Sex: M Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	08:39:26
Diagnostic Test	Pass	08:39:53
Air Blank	0.000	08:40:20
0.100 Cal Check	0.097	08:40:39
Air Blank	0.000	08:41:09
Subject Test	0.000	08:41:47
Air Blank	ABT*	08:41:53
Air Blank	0.000	08:42:22

\*Sequence Aborted

**Not a Successfully Completed Test Sequence**

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-000332  
04/01/2010 08:43:33

387  
On  
TS4

Standard Lot#: 642845  
Last Changed By: C. CONLEY #34517

QAS: C. CONLEY #34517  
TUCSON PD

Operator: C. CONLEY #34517  
TUCSON PD

Subject: TEST, SQAP  
DOB: 11/11/1111  
Sex: M Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	RFI*	08:44:05
Air Blank	IPS**	08:44:16

\*RFI Detect  
\*\*Improper Sample

**Not a Successfully Completed Test Sequence**

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-000332  
04/01/2010 08:45:13

457  
cu

Standard Lot#: 642845  
Last Changed By: C. CONLEY #34517

154

QAS: C. CONLEY #34517  
TUCSON PD

Operator: C. CONLEY #34517  
TUCSON PD

Subject: TEST, SQAP  
DOB: 11/11/1111  
Sex: M Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	08:45:58
Diagnostic Test	Pass	08:46:26
Air Blank	0.000	08:46:53
0.100 Cal Check	0.098	08:47:12
Air Blank	0.000	08:47:42
Subject Test	0.339*	08:50:50
Air Blank	0.000	08:51:24
Air Blank	0.000	08:51:51
Subject Test	XXX**	08:52:20
Air Blank	0.000	08:52:54

\*Deficient Sample

\*\*Invalid Sample

Begin new deprivation period

**Not a Successfully Completed Test Sequence**

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number: 80-000332  
04/01/2010 08:53:23

5067  
en

-----  
Standard Lot#: 642845  
Last Changed By: C. CONLEY #34517  
-----

TSA

-----  
QAS: C. CONLEY #34517  
TUCSON PD  
-----

Test	g/210L	Time
Air Blank	0.000	08:53:51
Cal Check	0.097	08:54:09
Air Blank	0.000	08:54:39

*Clare Conley*

Operator's Signature

EXHIBIT G-2

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

657  
Ch  
TH  
4.8.2010

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

QA Specialist Clare Conley AGENCY Tucson Police Department  
DATE 4-1-10 TIME 0837  
Intoxilyzer Serial # 80-000339 Location ODW DUI HQ

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.  
OR  
Pour a standard alcohol concentration solution \_\_\_\_\_ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 098 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Clare Conley

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number: 80-000332  
04/01/2010 08:37:00

7567  
Ln  
104

-----  
Standard Lot#: 642845  
Last Changed By: C. CONLEY #34517  
-----

QAS: C. CONLEY #34517  
TUCSON PD  
-----

Test	g/210L	Time
Air Blank	0.000	08:37:27
Cal Check	0.098	08:37:46
Air Blank	0.000	08:38:16

*Clare Conley*

\_\_\_\_\_  
Operator's Signature