



CERTIFICATION by CUSTODIAN of RECORDS

Tucson Police Department Crime Laboratory  
1306 W. Miracle Mile, Tucson, AZ 85705

TS4  
1-18-2012

I hereby certify that these documents are a true and correct copy of the periodic maintenance\* for the Intoxilyzer 8000 Serial Number 80-000334 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 1/5/12 Cal 1 0.098 AC Cal 2 0.097 AC

Between the dates of 12/8/11 and 1/5/12 the instrument was operating properly and accurately.

The instrument was taken out of service for repair on -.  
The instrument was returned to service on -.

Sarah Cione

Sarah L. Cione

1/11/12


ADPS QAS Permit # 33355

SUBSCRIBED AND SWORN to before me this 11, day of

January, 2012.

A. M. Hernandez

Notary Public

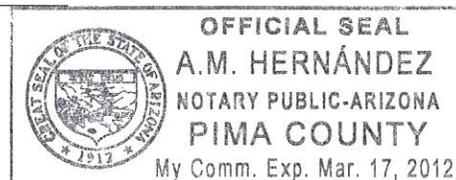

\*Includes documents specified in A.A.C. R13-10-104.A.5

1 of 7sc

EXHIBIT G-4

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

B4  
1-18-2012

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD QUALITY ASSURANCE PROCEDURE

QA SPECIALIST Sarah Cione AGENCY Tucson Police Department

DATE 1/5/12 TIME 1549

INTOXILYZER SERIAL # 80-000334 LOCATION DUI HQ

( ✓ ) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- ( ✓ ) 1. Clock time check.
- ( ✓ ) 2. Date check.

OPERATIONAL TESTS

- ( ✓ ) 1. Alcohol-free subject test result 0. 000 AC.
- ( ✓ ) 2. Error recognition logic system functioning.  
Not a Successfully Completed Test Sequence printed
- ( ✓ ) 3. Proper sample recognition system.  
Not a Successfully Completed Test Sequence printed  
Deficient sample printed.
- ( ✓ ) 4. Standard Calibration Check standard 0. 100 AC  
Result: 0. 097 AC

Instrument is operating properly and accurately. YES  NO

COMMENTS

---



---



---



---

SIGNATURE Sarah Cione

2 of 750

BSH

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-000334  
01/05/2012 15:50:39

Standard Lot#: 20910100A1  
Last Changed By: S. CIONE #50616

QAS: S. CIONE #50616  
TUCSON PD

Operator: S. CIONE #50616  
TUCSON PD

Subject: SQAP, SQAP  
DOB: 11/11/1111  
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	15:51:30
Diagnostic Test	Pass	15:51:58
Air Blank	0.000	15:52:25
0.100 Cal Check	0.099	15:52:44
Air Blank	0.000	15:53:13
Subject Test	0.000	15:53:37
Air Blank	0.000	15:54:04
Five Minute Wait	ABT*	
Air Blank	0.000	15:54:38

\*Sequence Aborted

### Not a Successfully Completed Test Sequence

3 of 7 sc

TS4

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-000334  
01/05/2012 15:55:34

Standard Lot#: 20910100A1  
Last Changed By: S. CIONE #50616

QAS: S. CIONE #50616  
TUCSON PD

Operator: S. CIONE #50616  
TUCSON PD

Subject: SQAP, SQAP  
DOB: 11/11/1111  
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	RFI*	15:55:56
Air Blank	IPS**	15:56:04

\*RFI Detect  
\*\*Improper Sample

### Not a Successfully Completed Test Sequence

4 of 75c

By

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number : 80-000334  
01/05/2012 15:56:31

Standard Lot#: 20910100A1  
Last Changed By: S. CIONE #50616

QAS: S. CIONE #50616  
TUCSON PD

Operator: S. CIONE #50616  
TUCSON PD

Subject: SQAP, SQAP  
DOB: 11/11/1111  
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

Test	g/210L	Time
Air Blank	0.000	15:57:12
Diagnostic Test	Pass	15:57:40
Air Blank	0.000	15:58:07
0.100 Cal Check	0.098	15:58:26
Air Blank	0.000	15:58:55
Subject Test	0.200*	16:02:04
Air Blank	0.000	16:02:37
Air Blank	0.000	16:03:05
Subject Test	XXX**	16:03:31
Air Blank	0.000	16:04:04

\*Deficient Sample  
\*\*Invalid Sample  
Begin new deprivation period

**Not a Successfully Completed Test Sequence**

5 of 7sc

BH

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number: 80-000334  
01/05/2012 16:08:15

-----  
Standard Lot#: 20910100A1  
Last Changed By: S. CIONE #50616  
-----

QAS: S. CIONE #50616  
TUCSON PD  
-----

Test	g/210L	Time
Air Blank	0.000	16:08:42
Cal Check	0.097	16:09:01
Air Blank	0.000	16:09:30

*Sarah Cione*

Operator's Signature

EXHIBIT G-2

6 of 7sc

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY  
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

STANDARD QUALITY ASSURANCE PROCEDURES  
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

TS4  
1.18.2012

QA Specialist Sarah Cione AGENCY Tucson Police Department

DATE 1/5/12 TIME 1549

Intoxilyzer Serial # 80-000334 Location DUIHQ

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.  
OR  
Pour a standard alcohol concentration solution \_\_\_\_\_ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 098 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Sarah Cione

DPS Form Exh G-2 (Rev. 05-01)

7 of 7 sc

INTOXILYZER 8000  
Location: TUCSON POLICE DEPT  
Serial Number: 80-000334  
01/05/2012 15:49:01

134

-----  
Standard Lot#: 20910100A1  
Last Changed By: S. CIONE #50616  
-----

QAS: S. CIONE #50616  
TUCSON PD  
-----

Test	g/210L	Time
Air Blank	0.000	15:49:28
Cal Check	0.098	15:49:47
Air Blank	0.000	15:50:17

*Sarah Cione*

Operator's Signature