



CS
9-27-10

CERTIFICATION by CUSTODIAN of RECORDS
Tucson Police Department Crime Laboratory
270 S. Stone Avenue Tucson, AZ 85701-1917

I hereby certify that these documents are a true and correct copy of the periodic maintenance* for the Intoxilyzer 8000 Serial Number 80-001163 maintained by the Tucson Police Department Crime Laboratory pursuant to the requirements of the Arizona Department of Public Safety for a quality assurance program.

A quality assurance specialist (QAS) tested the instrument on the following dates. The acceptable accuracy limits, as defined in A.A.C. R13-10-104.A.2, for a 0.100AC solution are 0.090AC to 0.110AC.

Date: 9.23.2010 Cal 1 0.098 AC Cal 2 0.099 AC

Between the dates of 8.26.2010 and 9.23.2010 the instrument was operating properly and accurately.

The instrument was taken out of service for repair on .
The instrument was returned to service on .

Terry Gallegos

Terry S. Gallegos


9-27-2010

ADPS QAS Permit # 33169

SUBSCRIBED AND SWORN to before me this 27, day of September, 20 10.

A.M. Hernandez

Notary Public

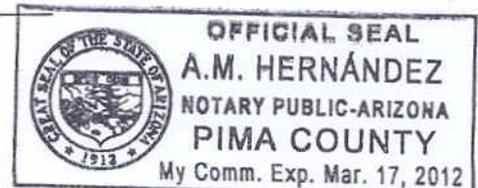

*Includes documents specified in A.A.C. R13-10-104.A.5

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY
STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

1057
TS4

STANDARD QUALITY ASSURANCE PROCEDURE

QA SPECIALIST Terry Gallegos 28847 AGENCY Tucson Police Department

DATE 9-23-2010 TIME 1400

INTOXILYZER SERIAL # 80-001163 LOCATION CrimeLab

(✓) 1. Display reads "PUSH BUTTON TO START"

DIAGNOSTIC TESTS

- (✓) 1. Clock time check.
- (✓) 2. Date check.

OPERATIONAL TESTS

- (✓) 1. Alcohol-free subject test result 0. 000 AC.
- (✓) 2. Error recognition logic system functioning. ✓✓✓
Not a Successfully Completed Test Sequence printed
- (✓) 3. Proper sample recognition system.
Not a Successfully Completed Test Sequence printed ✓
Deficient sample printed.
- (✓) 4. Standard Calibration Check standard 0. 100 AC
Result: 0. 099 AC

Instrument is operating properly and accurately. YES NO

COMMENTS

SIGNATURE

Terry Gallegos

CS

Zof 7
TS4

INTOXILYZER 8000
Location: ARIZONA
Serial Number : 80-001163
09/23/2010 14:02:17

Standard Lot#: 719919
Last Changed By: S. CIONE #50616

QAS: T. GALLEGOS #28847
TUCSON PD CRIME LAB

Operator: T. GALLEGOS #28847
TUCSON PD CRIME LAB

Subject: SQAP, SQAP
DOB: 11/11/1111
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 14:03:26 |
| Diagnostic Test | Pass | 14:03:53 |
| Air Blank | 0.000 | 14:04:20 |
| 0.100 Cal Check | 0.098 | 14:04:39 |
| Air Blank | 0.000 | 14:05:09 |
| Subject Test | 0.000 | 14:05:37 |
| Air Blank | ABT* | 14:05:47 |
| Air Blank | 0.000 | 14:06:15 |

*Sequence Aborted

Not a Successfully Completed Test Sequence

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

CS

3 of 7
B4

INTOXILYZER 8000
Location: ARIZONA
Serial Number : 80-001163
09/23/2010 14:06:22

Standard Lot#: 719919
Last Changed By: S. CIONE #50616

QAS: T. GALLEGOS #28847
TUCSON PD CRIME LAB

Operator: T. GALLEGOS #28847
TUCSON PD CRIME LAB

Subject: SQAP, SQAP
DOB: 11/11/1111
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 14:07:14 |
| Diagnostic Test | Pass | 14:07:41 |
| Air Blank | RFI* | 14:07:55 |
| Air Blank | IPS** | 14:08:06 |

*RFI Detect
**Improper Sample

Not a Successfully Completed Test Sequence

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

40 of 7
TS4

INTOXILYZER 8000
Location: ARIZONA
Serial Number : 80-001163
09/23/2010

14:08:20

Standard Lot#: 719919
Last Changed By: S. CIONE #50616

QAS: T. GALLEGOS #28847
TUCSON PD CRIME LAB

Operator: T. GALLEGOS #28847
TUCSON PD CRIME LAB

Subject: SQAP, SQAP
DOB: 11/11/1111
Sex: F Weight: 111

15 Minute Deprivation Period? Yes

| Test | g/210L | Time |
|-----------------|--------|----------|
| Air Blank | 0.000 | 14:09:13 |
| Diagnostic Test | Pass | 14:09:40 |
| Air Blank | 0.000 | 14:10:08 |
| 0.100 Cal Check | 0.098 | 14:10:26 |
| Air Blank | 0.000 | 14:10:56 |
| Subject Test | 0.252* | 14:14:04 |
| Air Blank | 0.000 | 14:14:37 |
| Air Blank | 0.000 | 14:15:04 |
| Subject Test | XXX** | 14:15:21 |
| Air Blank | 0.000 | 14:15:54 |

*Deficient Sample

**Invalid Sample

Begin new deprivation period

Not a Successfully Completed Test Sequence

18

5077
TS4

INTOXILYZER 8000
Location: ARIZONA
Serial Number: 80-001163
09/23/2010 14:16:13

Standard Lot#: 719919
Last Changed By: S. CIONE #50616

QAS: T. GALLEGOS #28847
TUCSON PD CRIME LAB

| Test | g/210L | Time |
|-----------|--------|----------|
| Air Blank | 0.000 | 14:16:40 |
| Cal Check | 0.099 | 14:16:59 |
| Air Blank | 0.000 | 14:17:29 |



Operator's Signature

THIS REPORT PREPARED PURSUANT TO DUTY IMPOSED BY
A.A.C. R13-10-104(A)

ARIZONA DEPARTMENT OF PUBLIC SAFETY

6087
TS4

STANDARD QUALITY ASSURANCE PROCEDURES
INTOXILYZER MODEL 8000

STANDARD CALIBRATION CHECK PROCEDURE

QA Specialist Terry Gallegos 28847 AGENCY Tucson Police Department
DATE 9-23-2010 TIME 1400
Intoxilyzer Serial # 80-001163 Location Crime Lab

- (✓) 1. Ensure that gas tank is attached to instrument and contains a standard alcohol concentration solution 0. 100 AC.
OR
Pour a standard alcohol concentration solution _____ AC, into a clean dry simulator and assemble the simulator. Ensure that a tight seal has been made. Turn on the simulator and allow temperature to reach 34°C ± 0.2°C
- (✓) 2. Intoxilyzer 8000 display reads "PUSH BUTTON TO START"
- (✓) 3. Go to the "Control Testing Menu". Select "D" for dry control test or "W" for wet control test. After selection is made press ENTER.
- (✓) 4. Air blank completed.
- (✓) 5. Calibration check completed. Test results 0. 098 AC.
- (✓) 6. Air blank completed.
- (✓) 7. Remove printed record. Attach the record to the completed checklist.

SIGNATURE Terry Gallegos

Pursuant to section 28-1327, Arizona Revised Statutes, this document is a certified duplicate of the information contained in computer storage devices of the Tucson Police Department.

CS

INTOXILYZER 8000
Location: ARIZONA
Serial Number: 80-001163
09/23/2010 14:00:39

7of7
156

Standard Lot#: 719919
Last Changed By: S. CIONE #50616

QAS: T. GALLEGOS #28847
TUCSON PD CRIME LAB

| Test | g/210L | Time |
|-----------|--------|----------|
| Air Blank | 0.000 | 14:01:07 |
| Cal Check | 0.098 | 14:01:25 |
| Air Blank | 0.000 | 14:01:55 |



Operator's Signature